


MODEL S15 Insert Earphones for fMRI Research Wiring Diagram

Male Banana Plug Or
Bare Wire Connection

Transformer

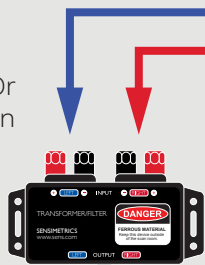


Patch Cables With
Male BNC 50Ω Plugs
On Each End
(Not Included)

Transformer Ground Cable

Audio Amplifier
(Not Included)



PC / Signal Source
(Not Included)

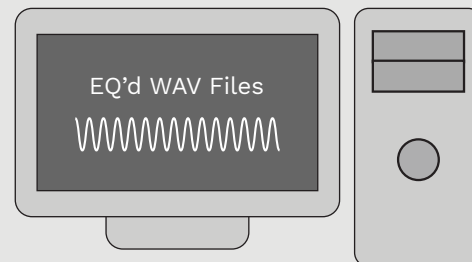

Penetration Panel

Direct Ground Connections
To Faraday Shield Ground

SCANNER ROOM

Shield Cable
(Available Separately)

Long Cable Assembly

~9 Meters

Filter

S15 Insert Earphones

~ 1.4 Meters

